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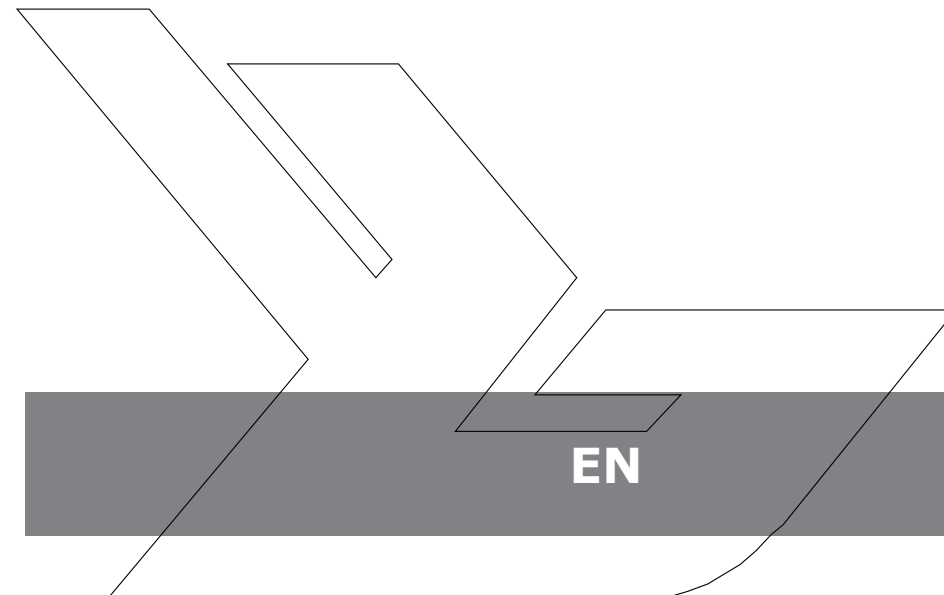
Sealed Lead Acid
Battery
Must be recycled
Properly

USER MANUAL
PRECISION SCALE

1. INTRODUCTION

Thank you for choosing our product. Each scale has a range of options and features that are ideal both for use in laboratory, classroom or for private use. We hope you enjoy working with our balance!

2. PACKAGE CONTENTS

- scale
- Weighing platform
- Power Cable
- Manual

3. BEFORE FIRST USE

INSERTING THE BATTERIES

Remove the battery box cover and insert the batteries (AA)

SCREEN AND KEYS

- ON / OFF 4 *
- batteries Level 2 *
- NULL (Zero) is displayed when the zero option is activated 1 *
- STABIL displayed when the result is stable 3 *
- HHHHH: the weight is too big, the scale cannot weigh.

PLACEMENT OF THE DEVICE AND PRECAUTIONARY MEASURES

To keep your balance in a good condition, we recommend you follow these tips:

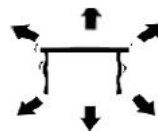
Avoid extreme temperatures. Do not place the unit in direct sunlight or near the air conditioning.

Make sure the scale is placed on a stable table and is not subject to vibrations.

Avoid unstable power sources. Do not use close to large electrical appliances such as welding machines or large motors. Do not leave batteries in the device if you do not use it for a long time.

Avoid vibrations. Do not place your scale near vibrating heavy machinery.

Avoid places where moisture can lead to condensation. Avoid direct contact with water.

Do not place the unit close to windows, doors or fans that may cause unstable results due to air currents.

Keep your machine clean. Do not store products on the scale when it is not used.

4. OPERATION

GENERAL INFORMATION

- a. Capacity / Size: 2000g / 0,01 g
- b. Display: LCD digital
- c. Backlight: blue
- d. Auto power save function: time off automatically programmable (2.5 or 8 min.)
- e. Units / Staging: g, ct, lb, oz, dr, GN, ozt, dwt, MM, tl.J, tl.T, tl.H, t (a total of 13 units)
- f. Counting function: 1. per unit 2. the scale can calculate a percentage
- g. Power: DC 9V adapter or AA batteries
- h. Weight: 550g

5. FUNCTIONS OF THE KEYS

ON / OFF

Press the ON / OFF button to start the unit. Hold the button pressed for 3 seconds to turn off the scale.

UNIT

The unit change is possible by pressing the UNIT button. Then press up to reach the desired unit.

PCS

Counting function: Press PCS to enter the counting function. The scales switches from the weighing mode to the counting mode. By pressing this button, you can switch from one mode to another.

ZERO

Press this button to subtract the weight of a container or any weight on the scale.

CALIBRATION AND PARAMETER SETTING

The calibration is performed in two stages. To calibrate the scale on 2kg, you need standards of 1 and 2 kg. Turn on the scale. Hold pressed the zero key for 3 seconds until the display shows „CALL“.

The identification number of the scale appears (SCL)

Press briefly again ON / OFF to display „CAL-1“. Wait until the scale displays „0“

Place the standard of 1 kg on the scale. It shows „PASS-1“ and „CAL-2“.

Then place the standard of 2 kg on the scale. It displays „PASS-2“

The scale is then calibrated.

FUNCTION OF TARE

1. Place an empty container on the weighing platform until the balance indicates a stable value. Then press „ZERO“ Your balance now displays „0“.
2. To exit the tare function, remove the container and repress the „ZERO“ button. Your scale is back to normal.

AUTO-ZERO

1. Press „ZERO“ if the scale does not show „0“ when nothing is placed on it.
2. The area of auto zero is $\pm 2\%$ of the total capacity of the scale.

DO NOT OVERLOAD THE BALANCE!

6. COUNTING MODE

1. Press „PCS“ to activate the counting function. Your weight unit is maintained.
2. Press „UNIT“, then choose between 100 and 1000 pieces and 10-1000 units by repeated pressure.
3. Place a unit / object on the scale. Press ON / OFF for the scale to remind the weight of a unit.

7. PERCENTAGE MODE

1. Press twice „PCS“ then press „UNIT“ to choose between 100.00 and 100.0
2. After choosing the percentage of accuracy, place the weight on the scale and press „ON / OFF.“ This weight now accounts for 100%. You can now place a weight, the scale displays the percentage.

NOTE

To return to weighing mode, press „PCS“

8. RESPECT THE WEEE (ENVIRONMENTAL STANDARD)

According to the European standard 2002/95/EEC, electrical or electronic components and must be recycled or disposed so that they do not damage the environment and the health. The recycling of batteries is clearly regulated by laws of 2002 and 2005 and was updated in the text WEEE (en: Recycling of electrical and electronic equipment.)

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expondo GmbH